# M291Ag-9 Silver Electrode

## Storage

Replace the protection cap, and store the electrode in its box at room temperature.

## M291Ag-9 Silver Electrode Electrode d'Argent

### Introduction

The M291Ag-9 Silver Electrode can be used directly or in a coated form for the determination of silver halides either by direct measurements or by potentiometric titration. It can be used together with a standard REF601 (Hg/Hg<sub>2</sub>SO<sub>4</sub>) reference electrode or with the REF251 or REF451 double junction reference electrode.

### Preparation for Measurement

1. Remove the protection cap from the electrode.
2. Before starting a measurement, rinse the electrode with distilled water. Take care not to wet the screw cap.
3. Check that the electrode body and sensing element are not damaged.

### Maintenance

#### Silver rod

The sensing part of the electrode can be polished with finishing abrasive bands, part no. B90X550 or silver polish.

#### Cleaning

The electrode must be carefully cleaned before use, but in many cases a thorough rinsing with distilled water will suffice. Owing to the high chemical resistance of the electrode, most cleaning chemicals and solvents may be used (the electrode tip and stem consist of pure silver and glass).

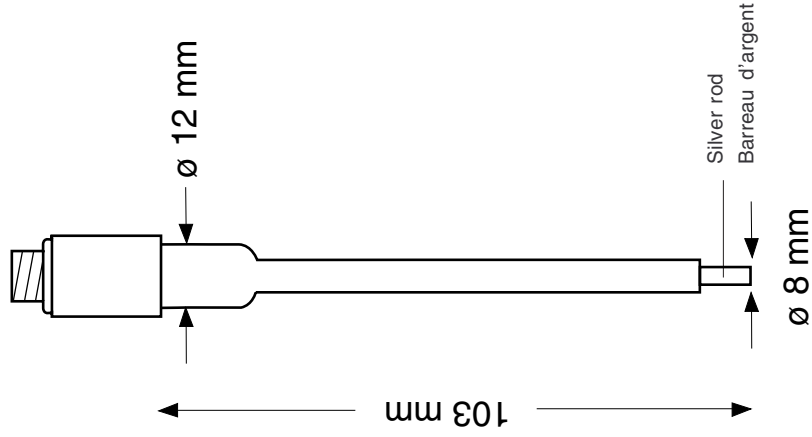
- To remove most electrode deposits which cannot be removed with pure water: flush with dilute nitric acid and/or ammonium hydroxide followed by distilled water.
- In case of severe contamination, clean the electrode with silver polish and then flush with acetone and distilled water.

### Accessories

BAO3 Finishing Abrasive Bands B90X550  
CL114 Electrode Cable, (1 m) with BNC plug A94L114  
CL116 Electrode Cable, (1 m) with type 7 plug A94L116  
CL111 Electrode Cable, (1 m) with banana plug A94L111

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D31M087 • Radiometer Analytical S.A. • France • 9809A
**RADIOMETER ANALYTICAL SAS**